

## Olympus microscope BX53-F



### To use Microscope

1. Place the sample on the bench below the lens of the microscope (keep the lens at a height while you do this).
2. Put on the lamp switch on the side of the microscope.
3. Watch the sample from the eyepiece.
4. Use focus, brightness, x and y position, lens magnification set knobs to obtain the required view.
5. Switch off the lamp.
6. Take off your sample.

### To take images

1. Put on the power switch on the panel (connecting the camera to the PC) on the table.
2. If the computer is switched off switch it on. Open the Analysis software if it is not already open.
4. Use the capture button and pull out the eye/camera rod (on the microscope) to capture the image you saw through the eyepiece.
5. Use the snapshot button to save the image on the workspace. Set magnification and take measurements.
6. Leave the software and computer on, switch off the power button on the panel on the table.
9. If you are the last user of the day, exit the software and shut down the computer.