

# Torrey-pines Low Temp Hot Plate

## Turn on procedure:

- Switch ON “Main switch”, Turn On backside switch of hotplate
- For setting up temp, Press “Temp button” then press “Down or Up” arrow key to decrease or increase temperature and again press “Temp button” to set temp.


Ex. Decrease Temp: Temp button → Down arrow key → Temp button.

Increase Temp: Temp button → Up arrow key → Temp button.

- For setting up time, Press “Timer button” then press “Down or Up” arrow key to decrease or increase time and again press “Timer button” to set time.

Ex. Decrease Time: Timer button → Down arrow key → Timer button.

Increase Temp: Timer button → Up arrow key → Timer button.

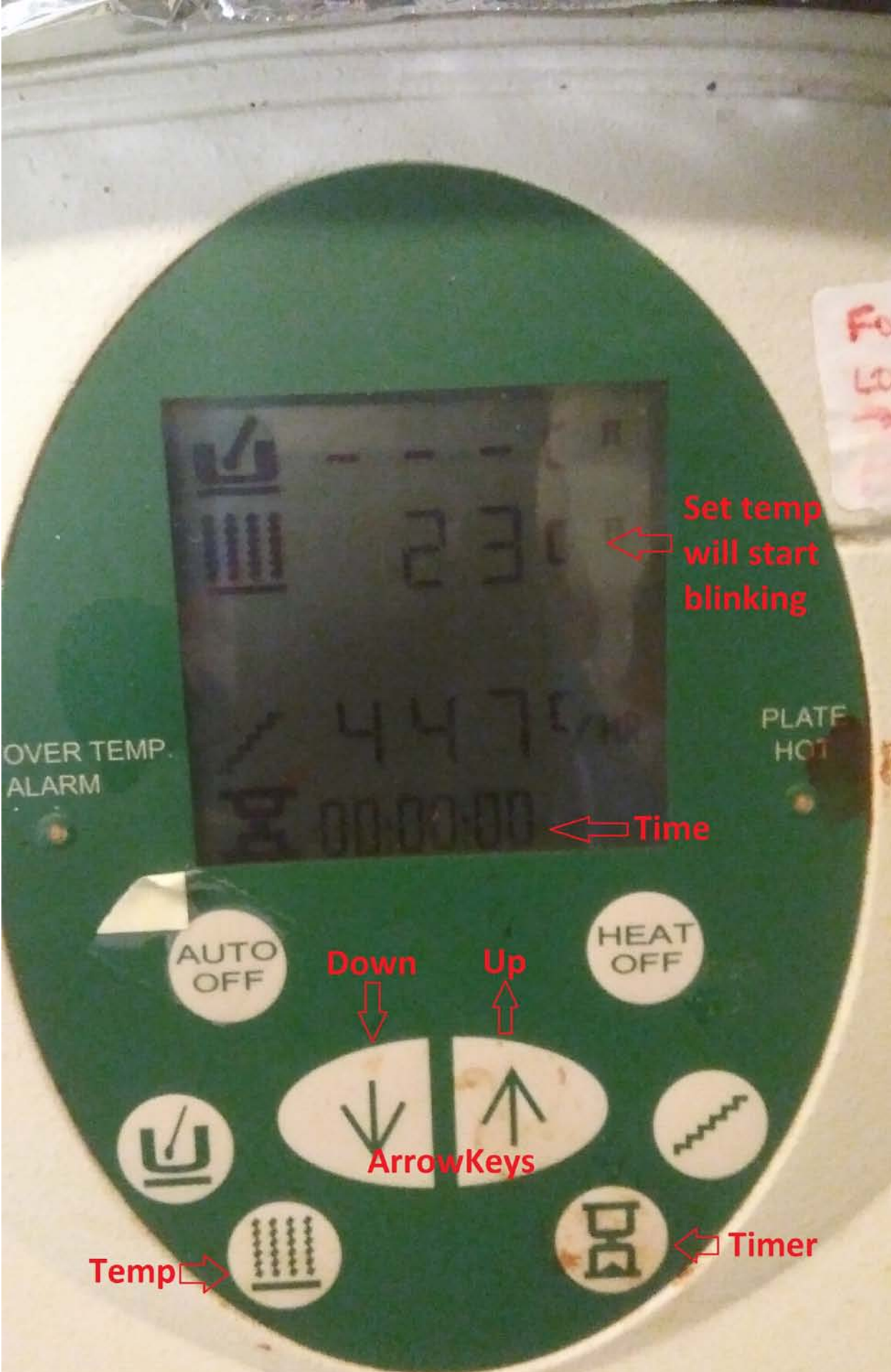

Set temp will start blinking

Time

Temp

Timer

Down

Up

ArrowKeys

OVER TEMP. ALARM

PLATE HOT

AUTO OFF

HEAT OFF

## **Turn OFF procedure:**

- Turn OFF backside switch of hotplate, Switch OFF “Main switch”.